

REDUCING END DETERMINATION THROUGH BICINCHONINIC ACID (BCA) ASSAY

Created/Updated by: Akanksha Nayak


Date: July/2014

Moran-Mirabal Research Group,
Department of Chemistry and Chemical Biology,
McMaster University
Hamilton, ON, Canada

<http://www.chemistry.mcmaster.ca/moran-mirabal>

BACKGROUND

- THE reducing end of a carbohydrate is a carbon atom that can be in equilibrium with the open-chain aldehyde or ketone form.

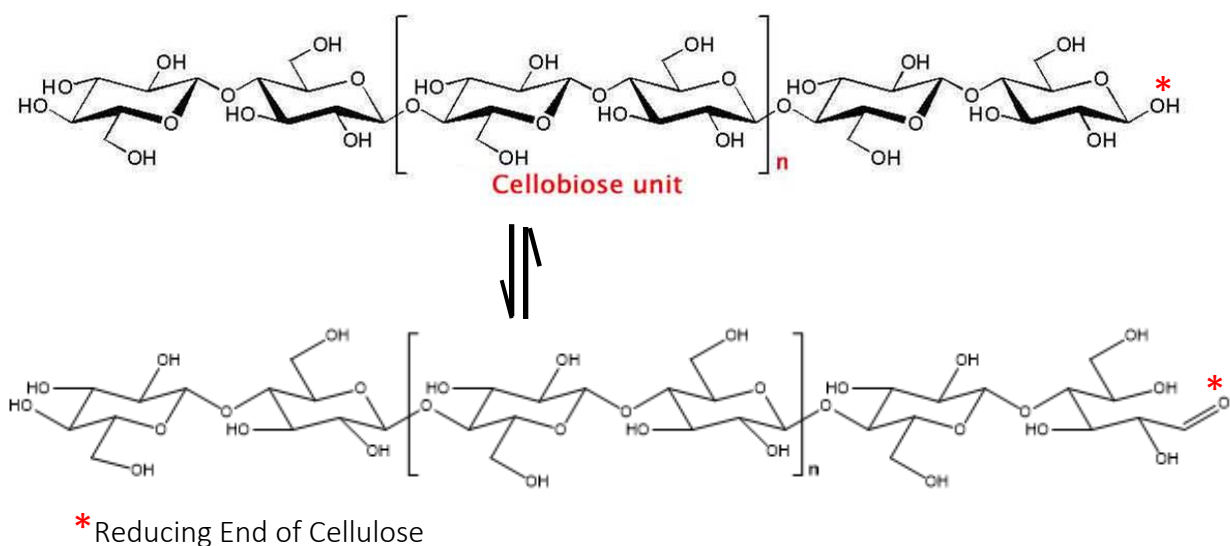

- This protocol determines the reducing end concentration of carbohydrate or cellulose samples.
- A temperature dependent reaction occurs where the reducing ends in the carbohydrate sample reduce Cu^{2+} ions from the copper(II) sulfate to Cu^+ . The amount of Cu^{2+} reduced is proportional to the amount of reducing ends present in the solution. Next, two molecules of bicinchoninic acid chelate with each Cu^+ ion, forming a lavender-colored product that strongly absorbs light at a wavelength of 560 nm.

A. MATERIALS

- BCA disodium salt hydrate
- Copper(II) sulfate pentahydrate ($\text{CuSO}_4 \cdot 5\text{H}_2\text{O}$)
- Glass Dish (pyrex)
- Graduated cylinder (pyrex or quartz)
- Greiner 96 well micro plate
- Hot plate with electronic stirring

REDUCING END DETERMINATION THROUGH BICINCHONINIC ACID (BCA) ASSAY

- L-serine
- Magnetic, Teflon coated stir bar
- Measuring balance
- Metal Stand
- Milli-Q
- Sodium bicarbonate (NaHCO_3)
- Sodium carbonate (Na_2CO_3)
- Sorvall Micro21R Centrifuge (Thermo Scientific)
- Styrofoam micro centrifuge stand
- Thermometer that goes up to 100°C
- Vortex Mixer (Fischer Scientific)
- 1.5mL Micro centrifuge tubes

B. PROTOCOL

1. Prepare Solution A with a pH of 9.7 by dissolving 27.14g of Na_2CO_3 , 12.1g of NaHCO_3 and 0.971g of BCA disodium salt hydrate in 500 mL distilled water.
2. Prepare Solution B with a pH of 3.4 by dissolving 0.624g of $\text{CuSO}_4 \cdot 5\text{H}_2\text{O}$ and 0.631g of L-serine in 500ml of distilled water.
3. Freshly prepare the BCA reagent by mixing equal volumes of solution A and solution B for each assay.
4. Mix equal amounts of the BCA reagent and sample. The amounts and concentrations of samples used are as follows:
 - BMCC- 0.5mL of 6 mg/mL (approx.)
 - Avicel- 0.5mL of 5 mg/mL
 - CF-11- 0.5mL of 5 mg/mL
 - CNC- 0.5mL of 2 mg/mL
5. After agitation using a vortex mixer, incubate the tubes at 75°C for 30 min in a water bath (prepared using the glass dish, heating plate and magnetic stirrer).
6. Cool the tubes for about an hour at room temperature.
7. Agitate the samples using a vortex mixer.
8. For insoluble samples, separate by centrifugation for 2 min at a speed of 10,000g.
9. Measure the absorbance at 560nm using a 96 well micro plate.

Note:

1. Solutions A and B are stable for at least a month at 4°C .
2. $1\mu\text{M}$ to $70\mu\text{M}$ glucose or cellobiose can be used as standards.


C. DISPOSAL

1. Pour the solution in the appropriate chemical waste container.

REDUCING END DETERMINATION THROUGH BICINCHONINIC ACID (BCA) ASSAY

D. SAMPLE RESULTS

Sample Calibration Curve with Reducing End Concentration of BMCC:

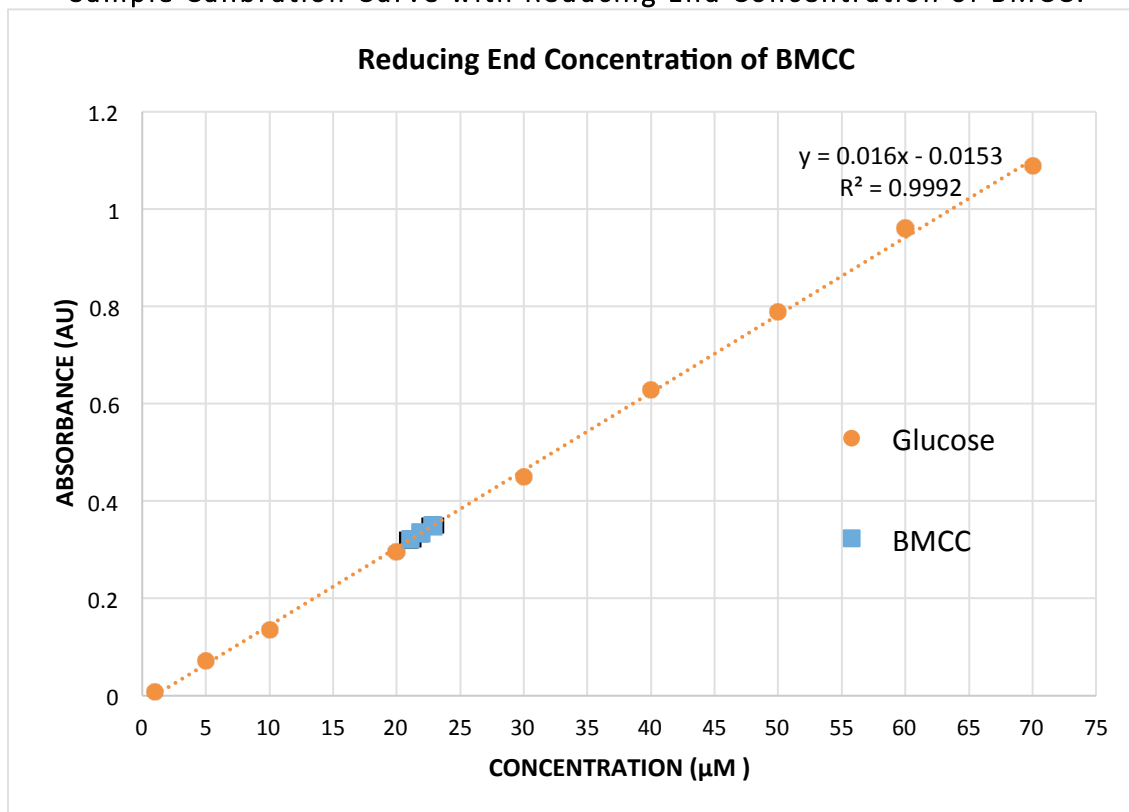

Examples of Reducing End Concentration Measured for Cellulosic Substrates:

Name	Reducing end (umol/g)	Error
Avicel PH-101	9.304377247	0.181488
	9.101460602	0.270536
	9.408505169	0.241984
BMCC (p-NDC)	3.165469349	0.111751
	3.429898844	0.115432
	3.286862141	0.081015
CNC (dil. 25%)	21.18618158	0.382342
	21.12062851	0.401058
	22.27209933	0.527542
CF-11	4.786769839	0.212963
	5.632591075	0.14403
	6.034831387	0.193498